SatPax® 70/30 IPA - Presaturated Wipes

70% IPA/30% WFI or DI Water

(en) Ready-to-use wipes for disinfection of bacteria and yeast on non-porous hard surfaces, floors, walls and ceilings in cleanrooms and other critical environments in pharmaceutical, biotechnology and healthcare industries, other life science applications and medical device manufacturing. Contains propan-2-ol (70% v/v): highly flammable liquid and vapour. Store in original container, protected from direct sunlight in a dry, cool and well-ventilated place, away from incompatible materials, food and drink; keep container tightly closed and sealed until ready for use. Do not freeze. **Directions for use:** Use at room temperature. Clean visibly soiled surfaces prior to disinfection. Remove wipe(s) from pack. Close Pack. Quarter fold the wipe (for mopping attach 4 layers on the mop head) and refold (or rotate) as needed to use each surface of the wipe(s). Apply wipe(s) directly to the surface to be disinfected. Use sufficient wipes to maintain a wet surface for a contact time of 1 minute for bacteria and yeast. After use immediately dispose of wipe(s) in a closed container. Professional use only. Refer to Summary of Product Characteristics (SPC) for additional information.

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